



Using Indicator Solutions

1. Add a few drops of each solution (acid, neutral or alkali) in the labelled column **(Do not mix the droppers!)**
2. To each dimple add a couple drops of indicator
3. Record the results

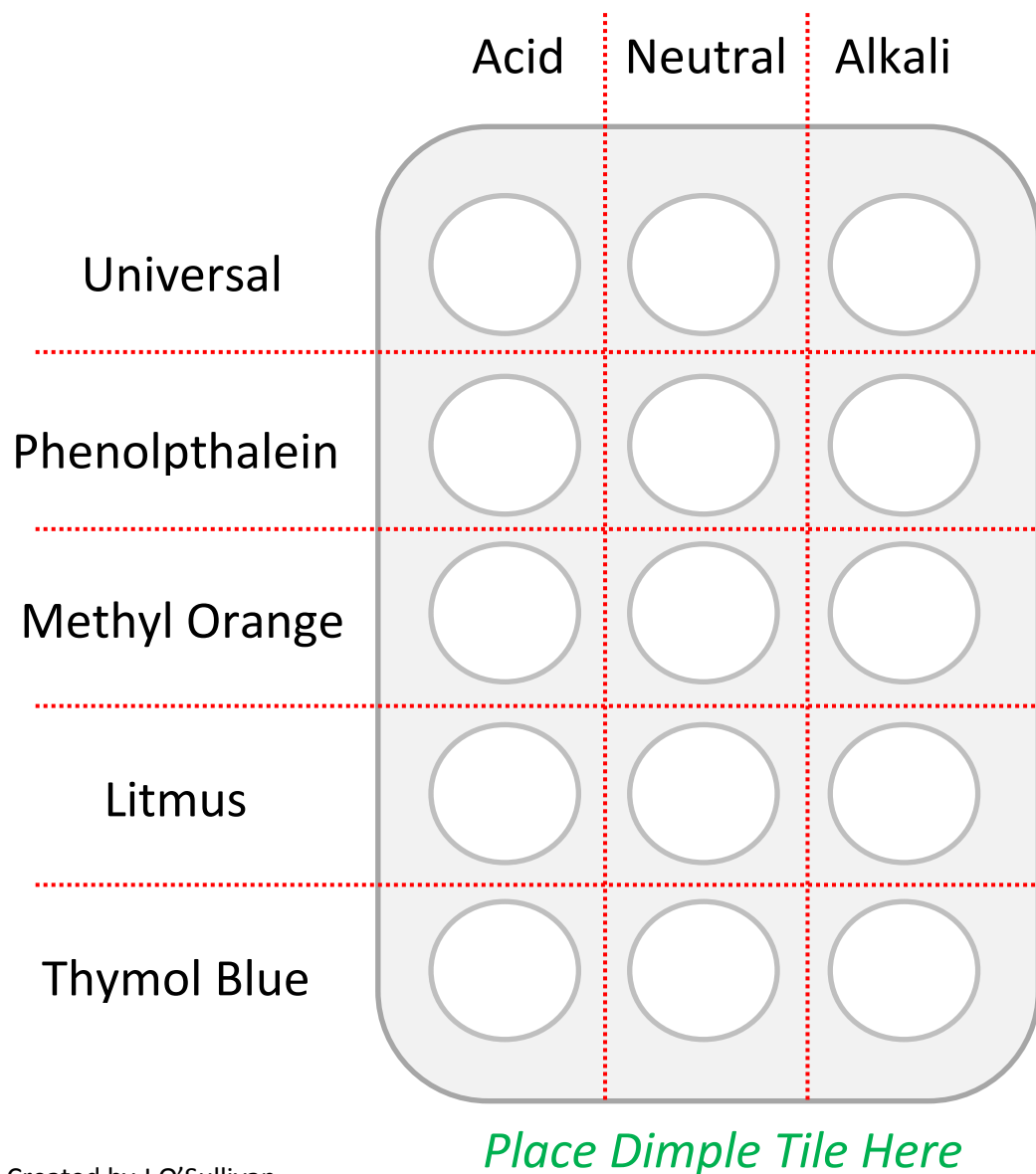

Using Indicator Solutions

1. Add a few drops of each solution (acid, neutral or alkali) in the labelled column **(Do not mix the droppers!)**
2. To each dimple add a couple drops of indicator
3. Record the results

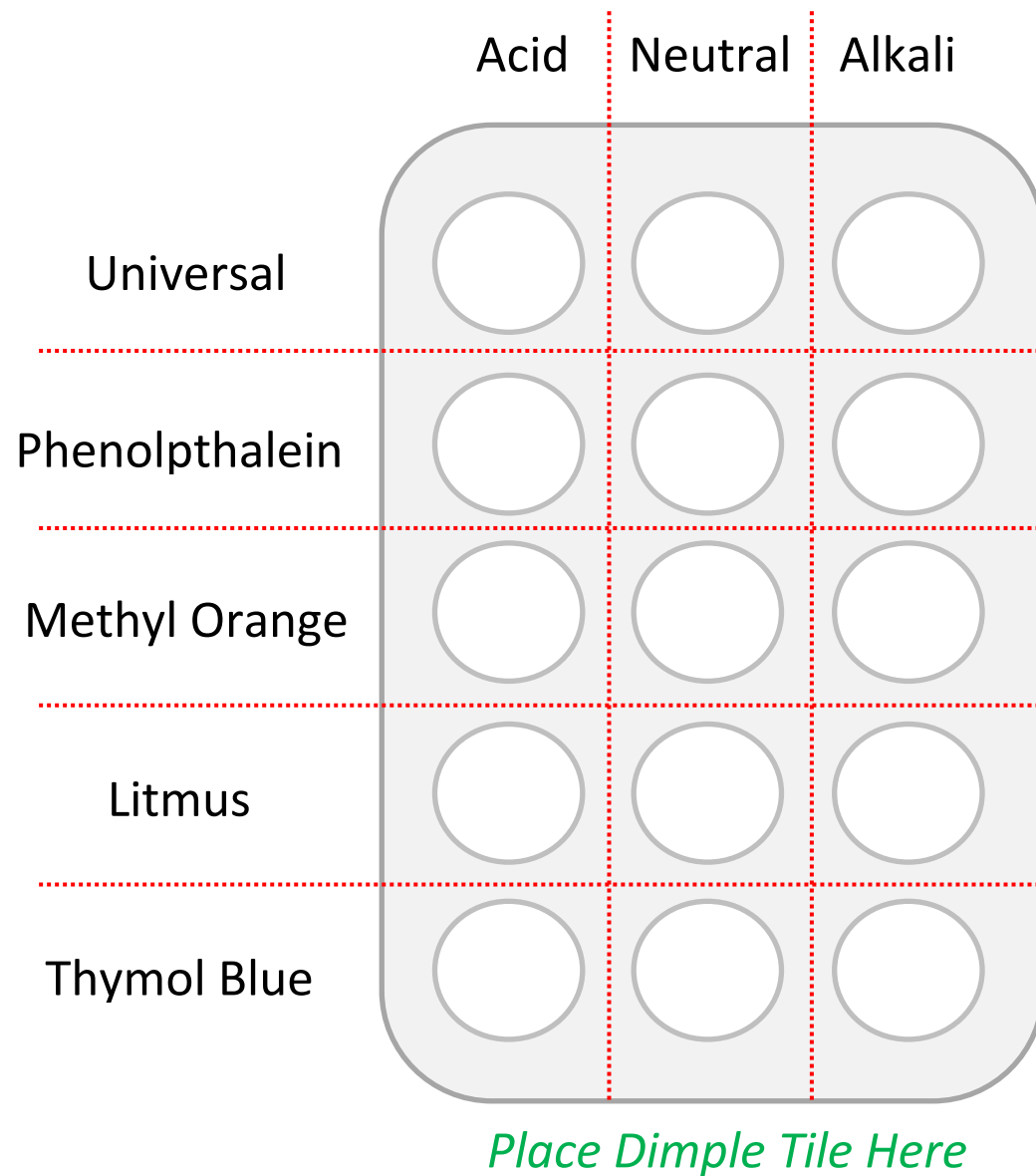

Indicator Solution	Colour		
	Acid	Neutral	Alkali
Universal			
Phenolphthalein			
Methyl Orange			
Litmus			
Thymol Blue			

Indicator Solution	Colour		
	Acid	Neutral	Alkali
Universal			
Phenolphthalein			
Methyl Orange			
Litmus			
Thymol Blue			

Indicator Solution	Colour		
	Acid	Neutral	Alkali
Universal			
Phenolphthalein			
Methyl Orange			
Litmus			
Thymol Blue			

Indicator Solution	Colour		
	Acid	Neutral	Alkali
Universal			
Phenolphthalein			
Methyl Orange			
Litmus			
Thymol Blue			

Indicator Solution	Colour		
	Acid	Neutral	Alkali
Universal			
Phenolphthalein			
Methyl Orange			
Litmus			
Thymol Blue			