



Step 1

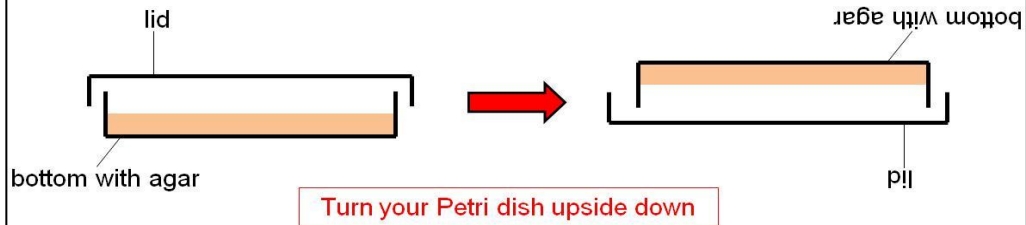


Step 2

A diagram of the inverted Petri dish with a red arrow pointing to the bottom surface, labeled 'bottom'. Below it, text reads: 'On the bottom of your Petri dish write your initials and the date'.

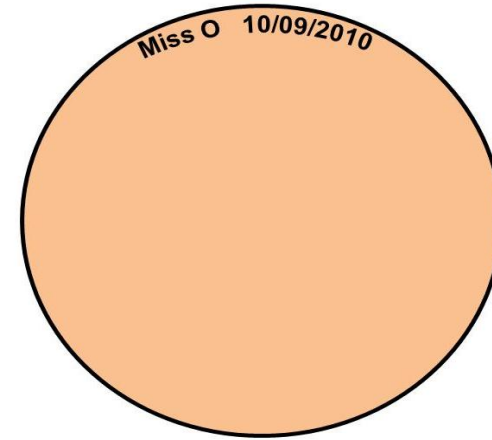


Step 3

A diagram of the inverted Petri dish with a red arrow pointing to the bottom surface, labeled 'bottom'. Below it, text reads: 'Mark the bottom of your Petri dish into quarters'.

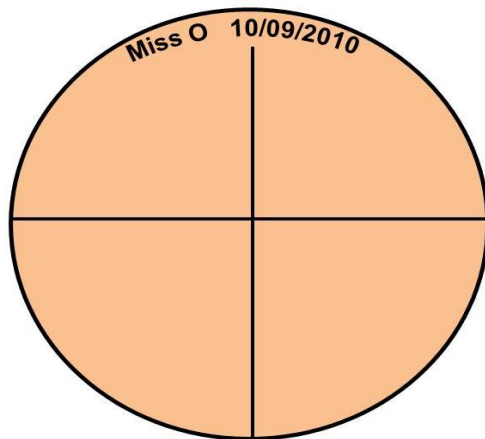


Step 4

A diagram of the inverted Petri dish with a red arrow pointing to the bottom surface, labeled 'bottom'. Below it, text reads: 'In each quarter write what you are swabbing'.

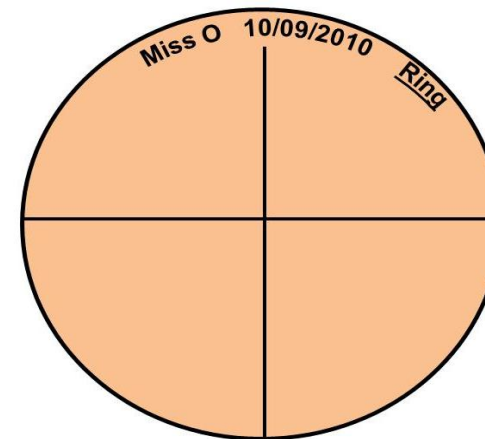

Step 5

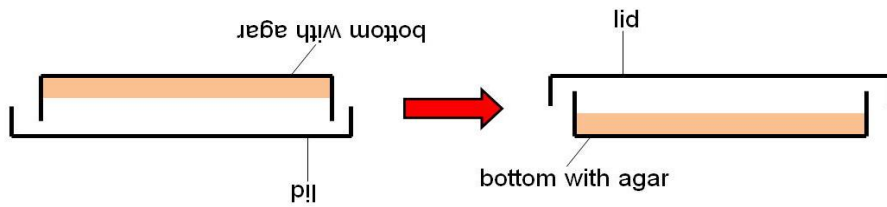


Flip over your Petri dish so the lid is back on top

Dip one of your cotton buds in water and rub what you want to test

Step 6

Take the lid off your Petri dish
Wiggle your cotton bud on the agar GENTLY

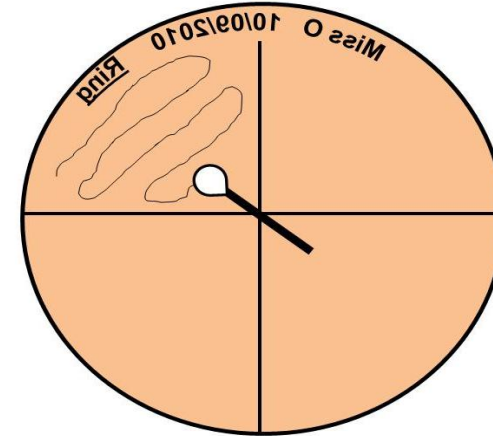


Remember to
put it in the
right quarter

Step 7




Put the lid back on your Petri dish between each test

Put the used cotton bud in Vircon and use a new one for the next test

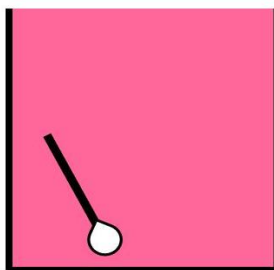

Step 8

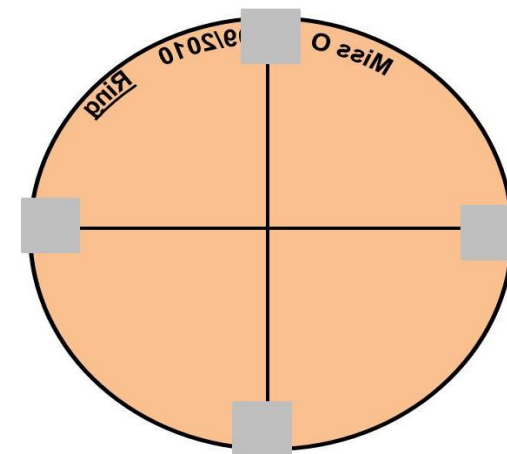

When you have tested 4 objects seal your dish with 4 pieces of sellotape